

Plasma-SeqSensei™ CRC RUO Kit



Master the sensitivity
in ctDNA detection

Plasma-SeqSensei™* (PSS) CRC RUO kits offer next-generation sequencing technology-based assays that enable highly sensitive and quantitative detection of mutations in circulating tumour DNA (ctDNA) from plasma and deliver results within two days as easy-to-read reports using Plasma-SeqSensei™ Software.

Key gene mutations covered include BRAF, KRAS, NRAS and PIK3CA to detect predictive markers, resistance mutations and frequently occurring clinically relevant genetic alterations in colorectal cancer research settings.

Unique benefits for clinical researchers



High sensitivity at low frequencies

PSS workflow reduces sequencing error rate up to 100-fold employing unique molecular identifiers (UID), enabling detection of 0.07% and higher mutant allele frequencies (MAF) with 95% certainty in a background of 10,000 wildtype copies.

→ Confident mutant variant reporting



Absolute mutant quantification

Internal quantifier Quantispik enables absolute ctDNA quantification down to a limit of detection (LOD) of seven mutant molecules (MM) independent of actual sample DNA input.

→ Consistent quantification independent of cfDNA input.

Unique benefits for clinical research laboratory



Short and standardised workflow

Standardised workflow from cfDNA samples to results in two days, including sequencing time.



Fast and convenient data analysis

Locally installed software automates data analysis and provides an easy-to-read mutation report, requiring only minimal bioinformatic knowledge.

Key facts

- ✓ Highly sensitive down to 0.07% MAF
- ✓ Able to detect six mutant molecules with 95% confidence
- ✓ Fast turnaround time (two days)
- ✓ Easy-to-read report



Target regions for PSS CRC RUO Kit

Gene ID#	Transcript ID#	CDS start	CDS end
BRAF	ENST00000288602	1,383	1,431
BRAF	ENST00000288602	1,742	1,813
KRAS	ENST00000256078	419	445
KRAS	ENST00000256078	169	228
KRAS	ENST00000256078	326	352
KRAS	ENST00000256078	34	102
NRAS	ENST00000369535	162	210
NRAS	ENST00000369535	420	449
NRAS	ENST00000369535	1	52
NRAS	ENST00000369535	341	364
PIK3CA	ENST00000263967	1,611	1,659
PIK3CA	ENST00000263967	3,118	3,195

Product specifications

Feature	Description
Starting sample	Whole blood and plasma
Sample capacity	2–16 samples per kit
QC function	Positive control and no template control (NTC) applied to every run
Input DNA required	5.7–95 ng / 116 µL
Number of amplicons	12
Sensitivity	0.07 % allele frequency with 95 % certainty in 10,000 wildtype copies
Cut-off	7 mutant molecules
Compatible sequencing instruments	Illumina MiSeq™ Illumina NextSeq 500/550™

* Plasma-SeqSensei™ CRC RUO Kit is for Research Use Only. Not for use in diagnostic procedures.